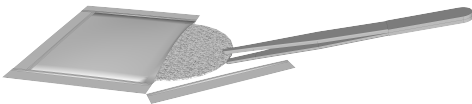
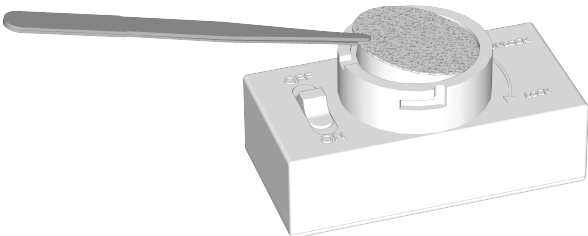
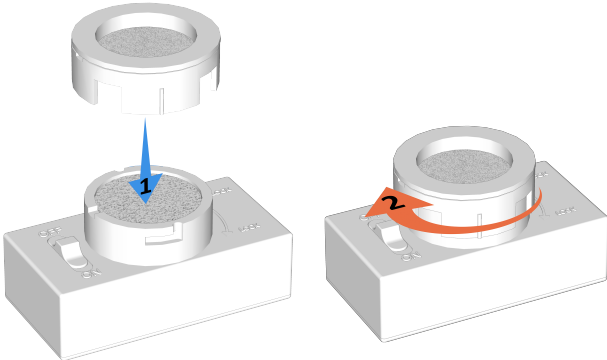
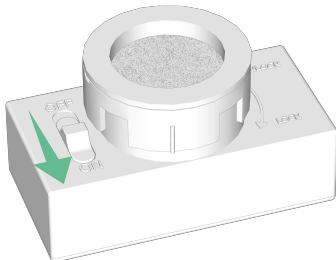


3	Open the package [6] and take the sample with the clamp [4]	
4	With the clamp [4], take a sampling sample [6] and insert it into the sampling device [2], positioning the pin in the slot provided for this purpose.	
5	Close the cover of the sampling device [2] by turning it clockwise until it clips	
6	Place the device where you want to make the sampling and switch the "ON / OFF" button to ON of the sampling device [2] and <b><i>Remember to note the time for choosing the duration of exposure</i></b>	
7	Toggle the "ON / OFF" button to OFF when the desired exposure time is reached. <b><i>Remember to note the end time of the exposure</i></b>	